Water distiller

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DESCRIPTION OF THE EQUIPMENT

CH. 2

- | | | |
|---------|------------|---|
| Pag. 32 | 2.1 | General description of the equipment |
| Pag. 32 | 2.2 | Table and description of the components |

TECHNICAL CHARACTERISTICS

CH. 3

- | | | |
|---------|------------|---|
| Pag. 32 | 3.1 | Technical characteristics and machine operating principle |
| Pag. 32 | 3.2 | Technical Data |

INSTALLATION AND FIRST-TIME USE

CH. 4

- | | | |
|---------|------------|-------------|
| Pag. 33 | 4.1 | Instruction |
|---------|------------|-------------|

MAINTENANCE

CH. 5

- | | | |
|---------|------------|----------|
| Pag. 34 | 5.1 | Cleaning |
|---------|------------|----------|

TROUBLESHOOTING

CH. 6

- | | | |
|---------|------------|------------------|
| Pag. 35 | 6.1 | Trouble shooting |
|---------|------------|------------------|

The data given in this manual may be modified without prior notice.

CHAPTER 1: GENERAL INFORMATION

1.1 TERMINOLOGY AND SYMBOLS

It is important that you understand the following symbols and their significance as they highlight important information such as critical situations, practical advice, or straightforward information.

If you have any doubts as to the significance of a symbol always consult this page.

- Critical situations concerning dangerous areas or tasks in which failure to observe the instructions may render the warranty null and void, or cause damage to equipment, connected components or the surrounding area, are highlighted by the symbol:

**WARNING!**

This symbol refers to safety standards which must be observed carefully in order to guarantee your safety, other people's safety and prevent damage to the equipment.

- Any general information and/or advice which may be useful at any time is highlighted by the following symbol:

**NOTE!**

This symbol highlights information and/or useful advice.

1.2 CONDITIONS OF GUARANTEE

This product is made of high quality material and great care has been taken in its manufacture. It is designed to give good service, provided it is properly operated and maintained. For your safety, please follow the instruction manual to prevent any possible damage.

The warranty shall not apply in the following circumstances: if it has been dismantled by an unauthorised person, connected to an incorrect mains voltage, or not used according to operating instructions, or to accidental damage.

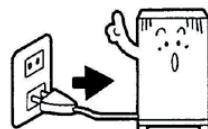
FOR YOUR SAFETY 1.3

Place the distiller on a dry, flat and heat-tolerant surface and keep it away from any heat source. Because the cooling fan uses air to cool it, we suggest you place it in a well-ventilated space.


Do not use a plastic bottle to store hot water.

Unplug the power line if you wish to terminate the distillation process prematurely.


Do not remove the upper cover when operating the distiller.

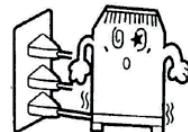Use a power point with proper voltage and do not share the same outlet with other electrical appliances to prevent overloading.

Do not let the children operate the distiller to avoid accidents.

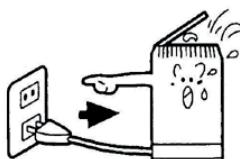Unplug the main power cord when filling or cleaning the distiller or when not in use.

Never immerse any part of the distiller in the water for the purpose of cleaning.

CHAPTER 2: DESCRIPTION OF THE EQUIPMENT

2.1 GENERAL DESCRIPTION OF THE EQUIPMENT

The water distiller makes the purest and cleanest water in the most natural and effective way.

Package includes:

Water distiller	Power cord
Protective cover	Residue cleaner
Activated charcoal (6 x)	Collection bottle

2.2 TABLE AND DESCRIPTION OF THE COMPONENTS

1. Upper cover
2. Cooling duct
3. Fan
4. Nozzle
5. Cushion rubber ring
6. Screws
7. Upper cover power line
8. Body
9. Power reset
10. Body socket

CHAPTER 3: TECHNICAL CHARACTERISTICS

3.1 TECHNICAL CHARACTERISTICS AND OPERATING START

When the distiller is turned on, the water temperature rises gradually ($\geq 100^{\circ}\text{C}$).

Chemicals, heavy metals and other contaminants, will be left as a residue at the bottom of the stainless steel tank. Impurities and unpleasant odours will be removed as steam collects on the surface of the cooling duct, which forms distilled water droplets and passes through the activated charcoal sachet. You then have sweet, pure water!

Please don't drink the first 2000cc.of water as this may contain any residues after manufacture.

3.2 TECHNICAL DATA

Power source	AC 100-240 / 50-60Hz
Power consumption	580 W (overall)
Capacity	4 lt
Fuse melting temperature	142 °C
Size	23 cm (Ø) x 36 cm (H)
Weight	3,5 Kg

CHAPTER 4: INSTALLATION AND FIRST-TIME USE

INSTRUCTION 4.1

Activated charcoal installation:

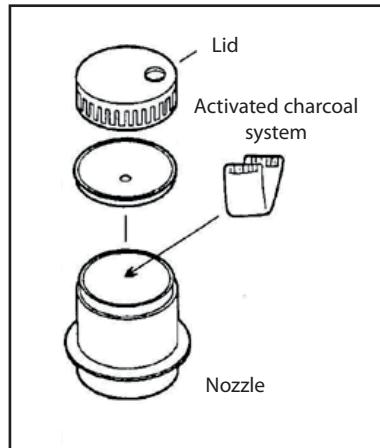
1. Turn the lid with the hole and separate it from the nozzle/square pouch.
2. Fold the activated charcoal sachet in half and put it into the nozzle and reassemble the lid.
3. Put the lid with the hole on to the collection bottle, lining it up with the spout on the distiller.

Each activated charcoal sachet can be used for up to a month.

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Pour water into the stainless steel tank

Use hot/cold water to fill the tank and make sure that you do not go beyond the full line. If you use hot water, you can shorten the time needed to distil water.

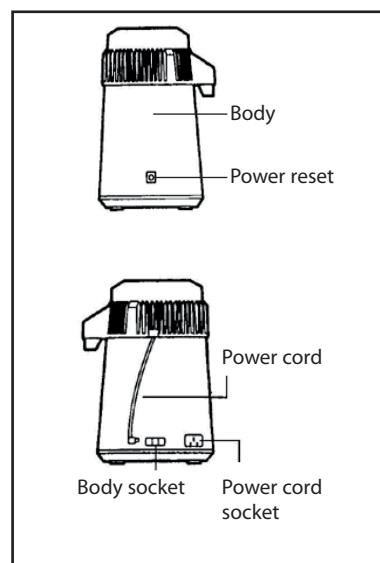
Starting up

Secure the upper cover and plug the power cord into the socket on the body. Place the collection bottle and the protective cover in a secure place and make sure they are levelled. Plug the power cord into the outlet and turn on the power switch (when first used the switch is automatically set in the "on" position, so there is no need to switch it on for the very first use).

You will start the process when turn the switch on and it will shut off automatically at the end of the process.

If you wish to terminate the process, unplug the power line from the outlet.

A power supply is turned off automatically when distillation is completed.

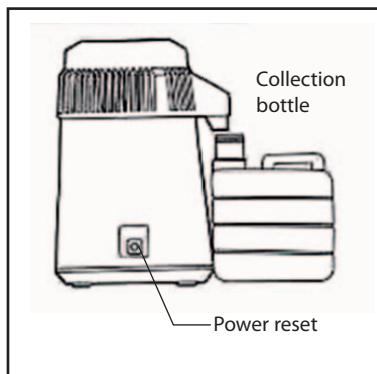
Approximately an hour after you have turned on the power switch, the distilled water begins to drip down into the collection bottle. It takes 5-6 hours to distil four litre of water.

When the power supply is cut off, you may move onto the next tank immediately.

Distilled water

For proper storage place the distilled water in the refrigerator in a sealed container. Before use shake the bottle to aerate the water.


**CHAPTER 5: MAINTENANCE****5.1 CLEANING**

Use household neutral detergent to clean the surface of the body and keep it dry (the rubber cushion within the upper cover is set with a specific direction; please do not remove it on your own).

To clean the body

The interior of the stainless steel tank is smooth. Never use abrasive cleaners on your distiller. After each cycle rinse the inside of the boiling chamber with clean water. To remove heavy deposits, use the distiller cleaner as follows:

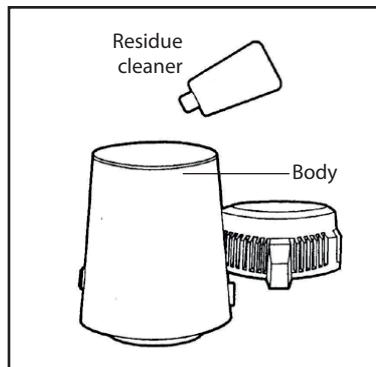1. Pour hot water into the stainless steel chamber until it covers the deposits.
2. Pour one to four spoonfuls of cleaner (based on the amount of deposits) into the tank.
3. Leave the upper cover off, plug in the main power plug, and turn on the power reset to heat the stainless steel tank until boiling.
4. Discard the water fifteen to twenty minutes later, then wash the interior with clean water.
5. Repeat the process until the chamber is clean.
6. Small amounts of residue are harmless, and you can continue to use it safely.

Cooling duct

If the condensing coil is clogged with dust, the efficiency of distillation may be affected. To clean the condensing coil, remove all three screws from the underside of the cover.

Use a tooth brush or a vacuum cleaner to remove the accumulated dust.

Others

"Take care" do not immerse the motor in the water when you are cleaning the cover. We recommend you regularly add a little oil on the motor shaft to help maintain the motor in good working order.

CHAPTER 6: TROUBLESHOOTING

5.1 TROUBLE SHOOTING

Conditions & Points to check:**Not making any distilled water:**

- Main power cord not plugged into power point or lid cord not plugged into base of the distiller.
- No power to electrical outlet.
- Reset button not pressed.

Not making enough distilled water during each cycle:

- Water not filled to maximum fill line.
- Lid gasket not sealing properly or up the wrong way.
- Collection bottle not placed directly under the spout.

Water or steam leaks from spout:

- Carbon sachets not properly placed in the nozzle.
- Collection bottle not correctly lined up with distiller.

Or contact your distributor for further advice.

This product must not be disposed of as household waste but rather, when no longer used, must be collected separately according to UE Directive 2012/19.

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